

NEBuilder[®] HiFi DNA Assembly

THE NEXT GENERATION OF DNA ASSEMBLY AND CLONING




be INSPIRED
drive DISCOVERY
stay GENUINE

Why choose NEBuilder HiFi

New to DNA Assembly?

NEBuilder HiFi DNA Assembly enables virtually error-free joining of DNA fragments, even those with 5' - and 3' -end mismatches. Available with and without competent *E. coli*, this flexible kit enables simple and fast seamless cloning utilizing a new proprietary high-fidelity polymerase. Find out why NEBuilder HiFi is the next generation of DNA assembly and cloning.

Not your average DNA assembly reagent

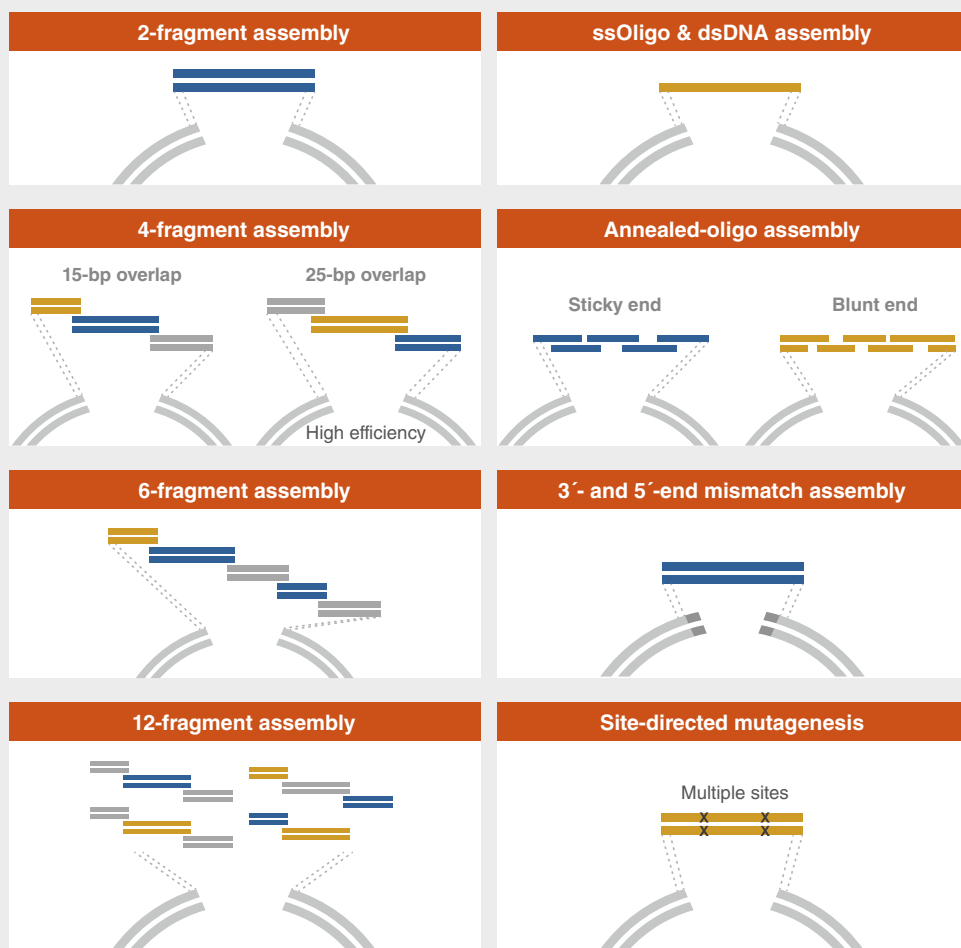

6 Reasons to choose NEBuilder HiFi

- 1 Save time**
Enjoy simple and fast seamless cloning in as little as 15 minutes.
- 2 Flexibility**
Use one system for both "standard-size" cloning and larger gene assembly products, up to 11 fragments.
- 3 Compatible with downstream applications**
DNA can be used immediately for transformation or as template for PCR or RCA.
- 4 Adaptable**
Adapts easily for multiple DNA manipulations, including site-directed mutagenesis.
- 5 Save money**
No licensing fees from NEB for NEBuilder products
- 6 Increased stability**
Store at -20°C, with improved stability over competition

Ordering Information


PRODUCT	NEB #	SIZE
NEBuilder HiFi DNA Assembly Master Mix*	E2621S/L/X	10/50/250 reactions
NEBuilder HiFi DNA Assembly Cloning Kit	E5520S	10 reactions
NEBuilder HiFi DNA Assembly Bundle for Large Fragments	E2623S	20 reactions

* Available in larger volumes. Please contact custom@neb.com for details.



DOWNLOAD THE NEB AR APP*

How does NEBuilder HiFi DNA Assembly work?

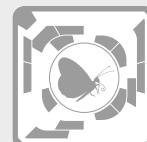

*see back cover for details

i DNA Assembly?


Using GeneArt® Gibson Assembly® or In-Fusion® Snap Assembly?

NEBuilder HiFi DNA Assembly offers several advantages over GeneArt Gibson Assembly and In-Fusion Snap Assembly. These include: higher accuracy due to the use of a high-fidelity polymerase, the ability to assemble both 5'- and 3'-end mismatches, lower DNA input requirements and the ability to bridge two double-stranded DNA fragments with a single-stranded DNA oligo (data not shown). NEBuilder HiFi DNA Assembly is the clear choice for efficient and accurate DNA assembly.

NEBuilder HiFi DNA Assembly offers improved efficiency in 4-fragment assembly reactions



4-fragment DNA assembly

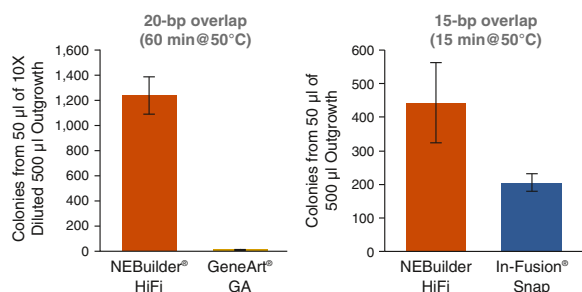


Four fragments (~20 fmol) with 20 bp overlap were assembled using NEBuilder HiFi DNA Assembly Master Mix (NEB #E2621), GeneArt Gibson Assembly Mix (Thermo Fisher® #A46627) and In-Fusion Snap Assembly Master Mix (Takara Bio USA #638947) to create a pUC19 vector. Assembly reactions were performed at 50°C for 60 min or 15 min. 2 µl of each assembled mix was transformed into NEB 5-alpha Competent E. coli (NEB #C2987) and spread on LB/Amp plates with IPTG and X-Gal. Blue colonies that indicated correct assembly were counted. NEBuilder HiFi DNA Assembly Master Mix yields more colonies than both competitors.

NEBuilder HiFi DNA Assembly offers improved efficiency in assembly of a ssDNA oligo with a linearized vector



ssOligo & dsDNA assembly

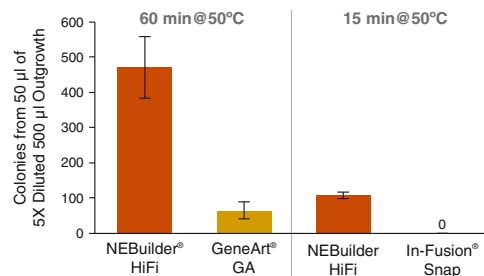

One pmol of ssDNA oligos were assembled with a linearized vector (30 ng of CRISPR Nuclease Reporter DNA) using NEBuilder HiFi DNA Assembly Master Mix (NEB #E2621), GeneArt Gibson Assembly Mix (Thermo Fisher #A46627) and In-Fusion Snap Assembly Master Mix (Takara Bio USA #638947). Assembly reactions were performed at 50°C for 60 min or 15 min. 2 µl of the assembled mix was transformed into NEB 5-alpha Competent E. coli (NEB #C2987). 20 colonies were further screened by PCR to confirm the presence of inserts. Greater than 95% of colonies tested from NEBuilder HiFi and GeneArt Gibson Assembly reactions contained proper inserts, although GeneArt Gibson Assembly yielded fewer colonies. In-Fusion Snap did not yield any successful colonies. NEBuilder HiFi DNA Assembly Master Mix outperformed both GeneArt Gibson Assembly and In-Fusion Snap Assembly.

Comparison of DNA Assembly Reaction Types

	NEBuilder HiFi DNA Assembly		GeneArt Gibson Assembly		In-Fusion Snap Assembly	
	Assembly efficiency	Covalently sealed?*	Assembly efficiency	Covalently sealed?*	Assembly efficiency	Covalently sealed?*
2-fragment assembly						
No mismatch	+++	Yes	++	Yes	++	No
3'- and 5'-end mismatch	+++	Yes	++	Yes	X	No
4-fragment assembly						
15-bp overlap & no mismatch	+++	Yes	++	Yes	++	No
25-bp overlap & no mismatch	+++	Yes	++	Yes	++	No
Oligo assembly						
3'- and 5'-overhang	+++	Yes	++	Yes	X	No
Blunt end & no mismatch	+++	Yes	++	Yes	X	No
ssOligo & vector	+++	Yes	NP	Yes	X	No

* Assembled products are treated with T5 exonuclease followed by PCR. Only covalently sealed products resistant to T5 exonuclease digestion can serve as templates for PCR and yield PCR product.

+++ Performs best; recommended
 ++ Performs well; but other product(s) perform better
 + Performs, but not recommended

X Does not perform
 NP Experiment not performed



Visit NEBuilderHiFi.com to view additional performance data, tutorials and much more



For help designing primers, try NEBuilder Assembly Tool at NEBuilder.neb.com

GERMANY & AUSTRIA
New England Biolabs GmbH
Brüningstr. 50, Geb B852
65926 Frankfurt/Main, Germany

Tel: +49/(0)69/305-23140
Fax: +49/(0)69/305-23149

Free Call: 0800/246 5227 (Germany)
Free Call: 00800/246 52277 (Austria)
info.de@neb.com

www.neb-online.de

HEADQUARTERS:

USA
New England Biolabs, Inc.

Telephone: (978) 927-5054
Toll Free (USA Orders): 1-800-632-5227
Toll Free (USA Tech): 1-800-632-7799

Fax: (978) 921-1350
info@neb.com
www.neb.com

www.neb.com



One or more of these products are covered by patents, trademarks and/or copyrights owned or controlled by New England Biolabs, Inc. For more information, please email us at busdev@neb.com. The use of these products may require you to obtain additional third party intellectual property rights for certain applications.

Your purchase, acceptance, and/or payment of and for NEB's products is pursuant to NEB's Terms of Sale at www.neb.com/support/terms-of-sale. NEB does not agree to and is not bound by any other terms or conditions, unless those terms and conditions have been expressly agreed to in writing by a duly authorized officer of NEB.

IN-FUSION® is a registered trademark of Takara Bio USA, Inc. GENEART® is a registered trademark of Thermo Fisher Scientific GeneArt GmbH LLC. THERMO FISHER SCIENTIFIC® is a registered trademark of Thermo Fisher Scientific, Inc. GIBSON ASSEMBLY® is a registered trademark of CODEX DNA, Inc. IPHONE® and IPAD® are registered trademarks of Apple, Inc.

Android™ is a trademark of Google, Inc.

© Copyright 2021, New England Biolabs, Inc.; all rights reserved.

Featured Tools



For help with finding the right products and protocols for each step of your next traditional cloning experiment try **NEBcloner®** at NEBcloner.neb.com



For help with designing primers for DNA assembly, try **NEBuilder® DNA Assembly Tool** at NEBuilder.neb.com



Download the **NEB AR App** for iPhone®, iPad® or Android™. Scan the augmented reality butterfly icon located on the corner of the page to find videos, tutorials and immersive experiences.

Your local NEB distributor:

BeNeLux:
BIOKE
Tel: (+31) 71 720 0220
(BE: 0800 - 71640)
info@bioke.com
www.bioke.com

BULGARIA
ELTA 90M Ltd.
Tel: (+359) 2 983 96 49
Fax: (+359) 2983 2211, 9274 232
office@elta90.eu
www.elta90.com

CROATIA:
Diagnostica Skalpeli d.o.o.
Tel: (+385) 1 377 8484
Fax: (+385) 1 377 8585
skalpeli@skalpeli.hr
www.skalpeli.hr

CZECH REPUBLIC:
BioTech a.s.
Tel: (+420) 2727 017 39
info@ibiotech.cz
www.ibiotech.cz

DENMARK:
BioNordika Denmark A/S
Tel: (+45) 39 56 20 00
Fax: (+45) 39 56 19 42
info@bionordika.dk
www.bionordika.dk

FINLAND:
BioNordika Oy
Tel: (+358) 207 410 270
info@bionordika.fi
www.bionordika.fi

GREECE:
BIOLINE SCIENTIFIC
Tel: (+30) 210 522 6547
Fax: (+30) 210 524 4744
info@bioline.gr
www.bioline.gr

HUNGARY:
Kvalitex Kft.
Tel: (+361) 340 4700
Fax: (+361) 339 8274
info@kvalitex.hu
www.kvalitex.hu

ITALY:
EUROCLONE S.P.A.
Tel: (+39) 02 38 1951
info@euroclone.it
www.euroclonegroup.it

ISRAEL:
Ornat BioChemicals &
Lab Equipment
Tel: (+972) 8 947 7077
Fax: (+972) 8 936 3034
ornatbio@ornat.co.il
www.ornat.co.il

LITHUANIA:
Nanodiagnostika, Ltd.
Tel: (+370) 525 052 44
lina@nanodiagnostika.lt
www.nanodiagnostika.lt

NORWAY:
BioNordika AS
Tel: (+47) 2303 5800
info@bionordika.no
www.bionordika.no

POLAND:
Lab-JOT Ltd. Sp.z o.o. Sp.k.
Tel: (+48) 2233 598 84
Fax: (+48) 2233 598 19
biuro@labjot.com
www.labjot.com

PORTUGAL:
Werfen
Tel: (+351) 214 247 300
dac-pt@werfen.com
www.pt.werfen.com

ROMANIA:
BioZyme
Tel: (+40) 264 523 281
contact@biozyme.ro
www.biozyme.ro

RUSSIA and CIS:
SkyGen LLC
Tel: (+88) 0033 312 26
info@skygen.com
www.skygen.com

SERBIA:
Alfa Genetics d.o.o.
Tel: (+38) 1011 2084 252
office@alfagenetics.rs
www.alfagenetics.rs

SPAIN:
Werfen
Tel: (+34) 900 822 290
customerservice-es@werfen.com
www.es.werfen.com

SLOVAK REPUBLIC
BioTech s.r.o.
Tel: (+421) 254 7744 88
Fax: (+421) 232 2010 48
info@ibiotech.sk
www.ibiotech.sk

SLOVENIA:
Mikro+Polo d.o.o.
Tel: (+386) 2614 33 00
Fax: (+386) 2614 33 20
podpora@mikro-polo.si
www.mikro-polo.si

SOUTH AFRICA:
Inqaba Biotechnical Industries
(Pty) Ltd
Tel: (+27) 1234 358 29
info@inqababiotech.co.za
www.inqababiotech.co.za

SWEDEN, ESTONIA, LATVIA,
LITHUANIA:
BIONORDIKA SWEDEN AB
Tel: (+46) 830 60 10
info@bionordika.se
www.bionordika.se

SWITZERLAND:
BioConcept Ltd.
Tel: (+41) 6148 680 80
Fax: (+41) 6148 680 00
info@bioconcept.ch
www.bioconcept.ch

TURKEY:
EKA BIOLAB TEKNOLOJİ
Tel: (+90) 312 248 13 24
Email: info@ekabiolab.com
www.ekabiolab.com

Distributors

For a complete list of all NEB subsidiaries
and distributors, please go to:

www.neb.com/international-support